


Reimagining DNA Payloads: Enzymatic Circular DNA for the Next Generation of Cell and Gene Therapies

Peter Archibald Ph.D, Mitchell Brun Ph.D., Elena Stoyanova Ph.D, Elena Garitta Ph.D, Duncan Simpson, Shivam Bhatt Ph.D, Jonathan Extance Ph.D, Tom Adie Ph.D, Milica Vukovic Ph.D MBA
Touchlight Genetics Ltd, Morelands & Riverdale Buildings, Lower Sunbury Road, Hampton, TW12 2ER



mbDNA demonstrates editing by both by EP and LNP

ElevateBio Authors: Aaron Hoover, Sharlene Amador, Chesney Michels, Jeff Cram, Mike Paglia

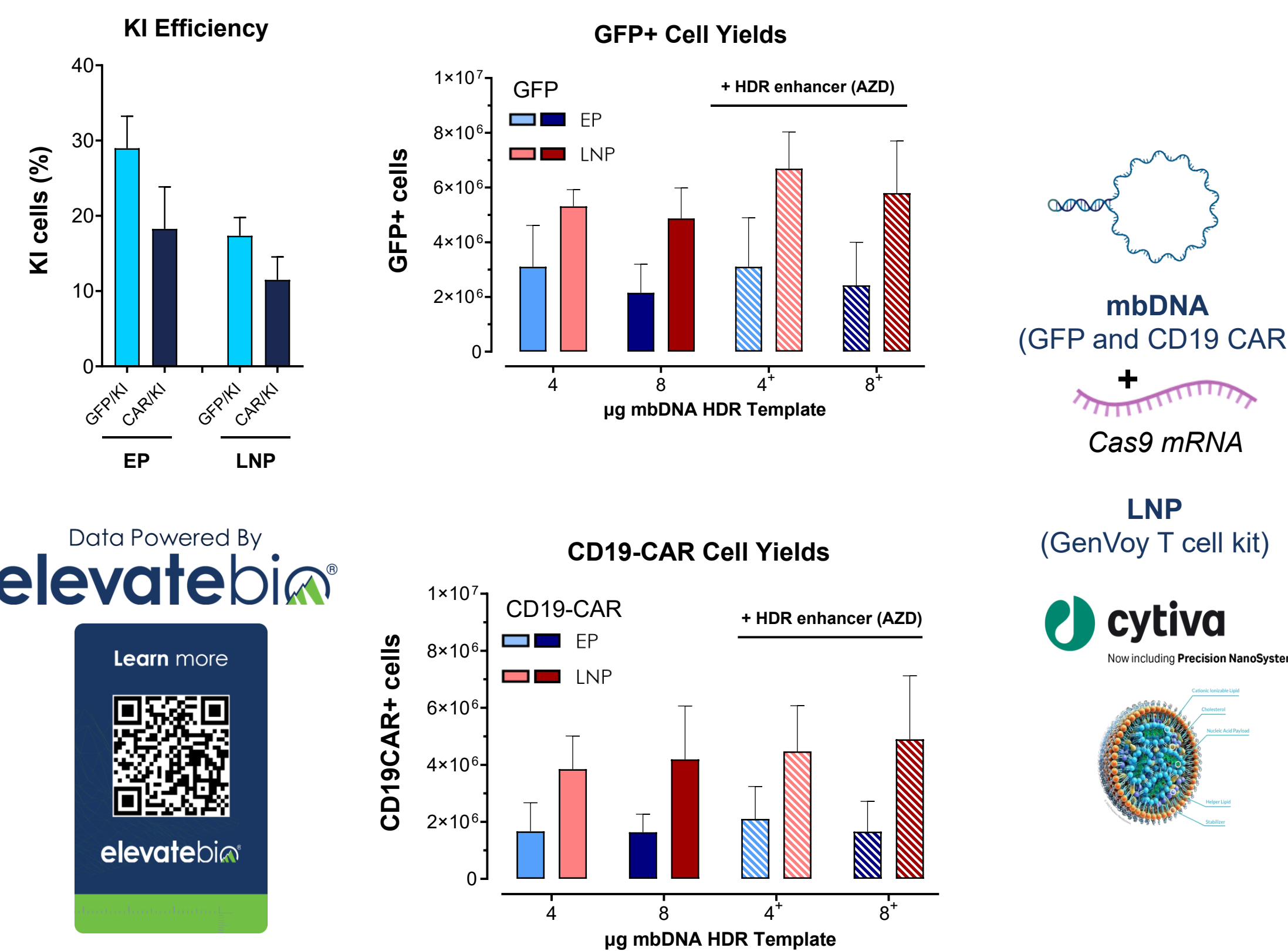


The superior expansion of **Cytiva LNP-treated cells** resulted in 1.7 to 3.0-fold higher total edited cell yields compared to EP-treated cells

LNP-formulated mbDNA+Cas9 mRNA + sgRNA successfully knock-in two different payloads in T cell ex vivo and achieve 17% GFP KI and 12% CAR KI

It is possible that higher KI efficiency can be achieved through process optimization

Improvement over next-gen pDNA with LEG14 nuclease

ElevateBio Authors: Aaron Hoover, Sharlene Amador, Chesney Michels, Jeff Cram, Mike Paglia

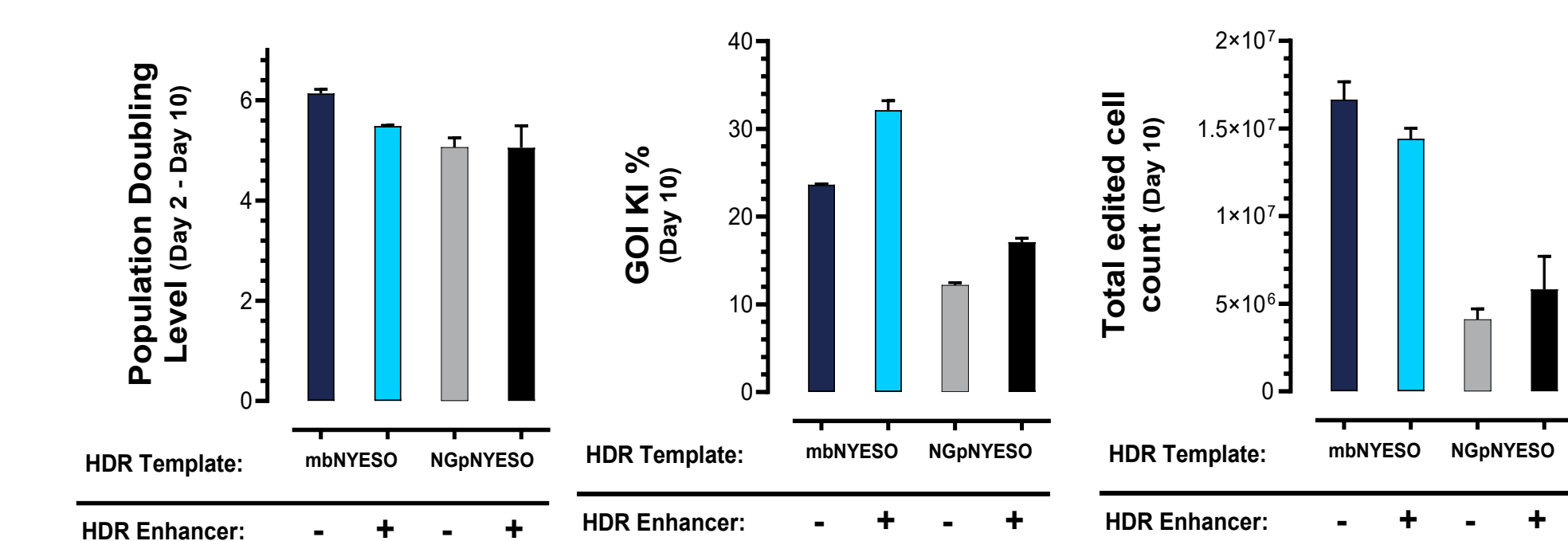


mbDNA outperforms NGpDNA with improved targeted TCR insertion of two donor templates with LEG14 nuclease

With ~4.5kb template, mbDNA achieved a 1.8-fold increase in HDR

With ~6.5kb template, mbDNA achieved a 2.8-fold increase in HDR


Data Powered By
elevatebio

mbDNA is tolerated well and has high KI

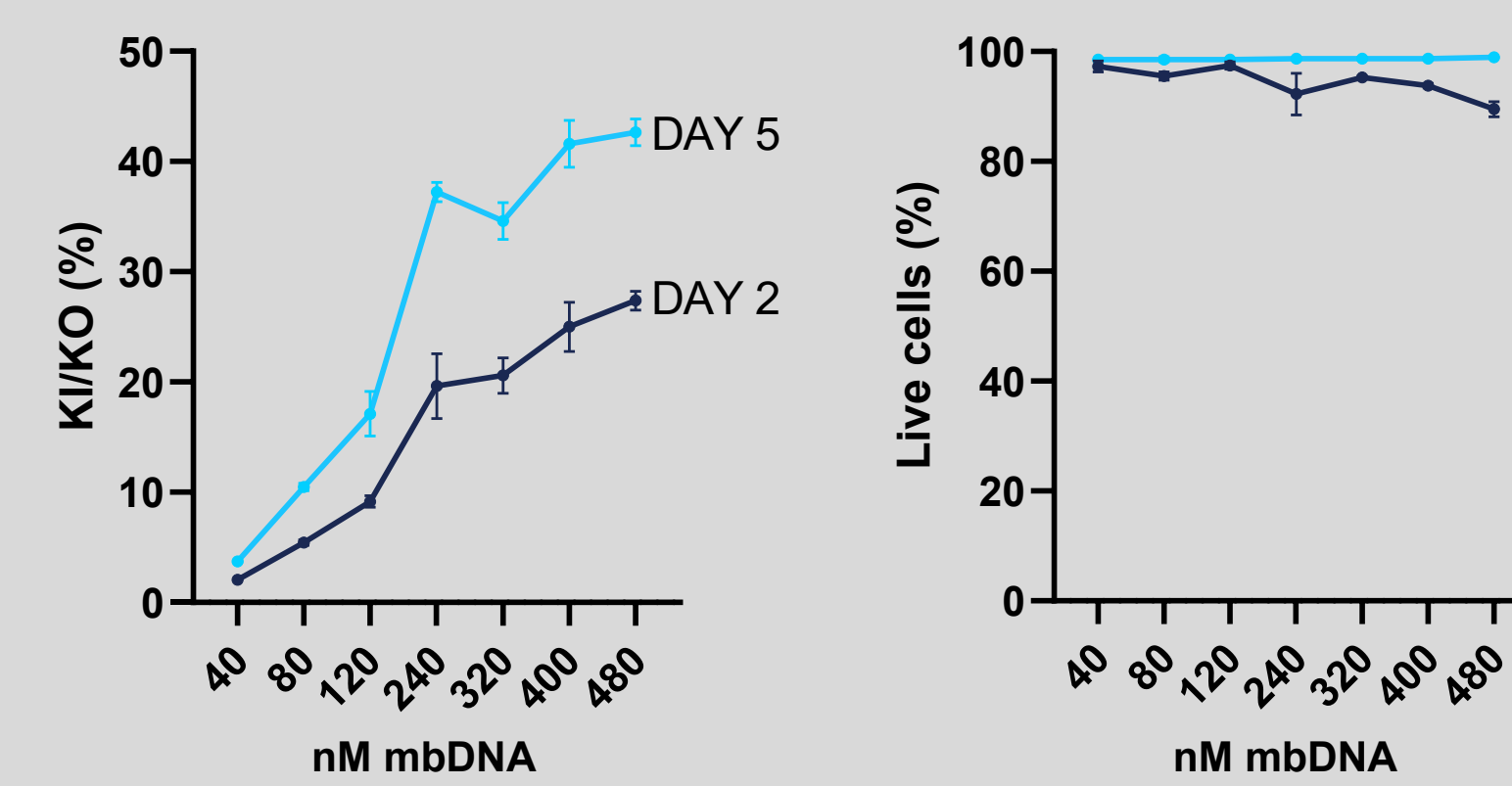

Toxic dose limit not found

Low donor-to-donor variability

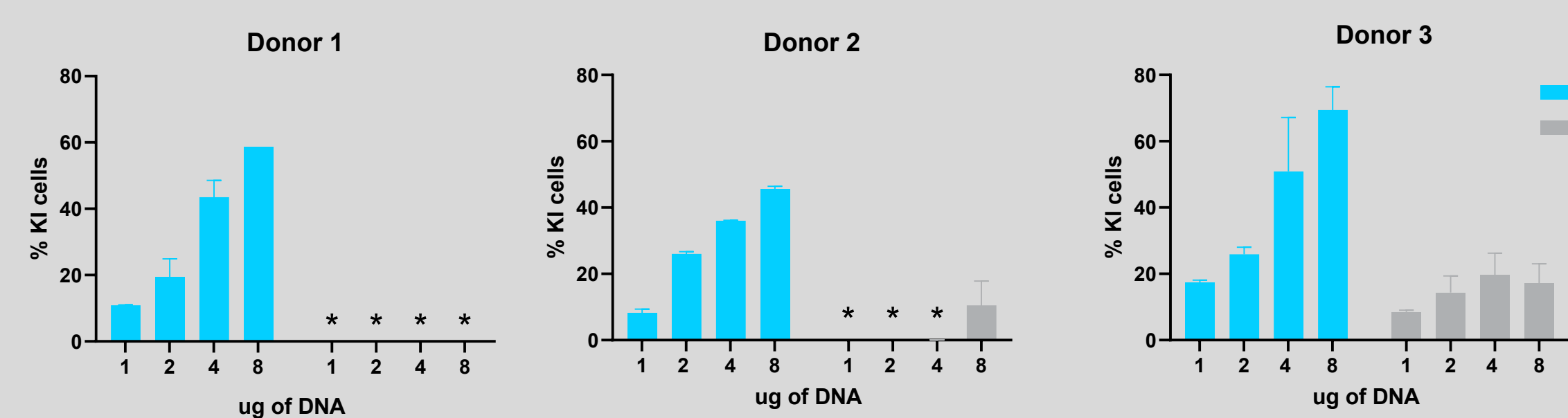

Consistent knock in across 3 donors

Multiple scales and batches performed similarly

Highly efficient editing produces CAR T cells

- 60-75% CAR KI achieved

Electroporation did not impair T cell expansion

- mbDNA samples expanded similarly to mock controls

High number of CAR⁺ T cells generated

- >1E9 CAR⁺ cells produced

Expected CD4:CD8 ratio maintained at all doses

- Low donor to donor variability observed

Low levels of terminally differentiated T cells were quantified post-electroporation

- CAR⁺CD57⁺% was low across conditions

T cells retain capacity for proliferation and survival post-electroporation

- CAR⁺CD8⁺CCR7⁺% maintained across conditions

- Synthetic **circular** ssDNA:
- Improved stability
- Structural advantage


- Gene-length:

- Multi-kb** sequences
- Beyond viral** packaging capacity

- Improved **manufacturability & scalability** over existing long ssDNA


- Leveraging Touchlight's established **enzymatic technology** and GMP manufacturing capabilities

Outperforms ssDNA

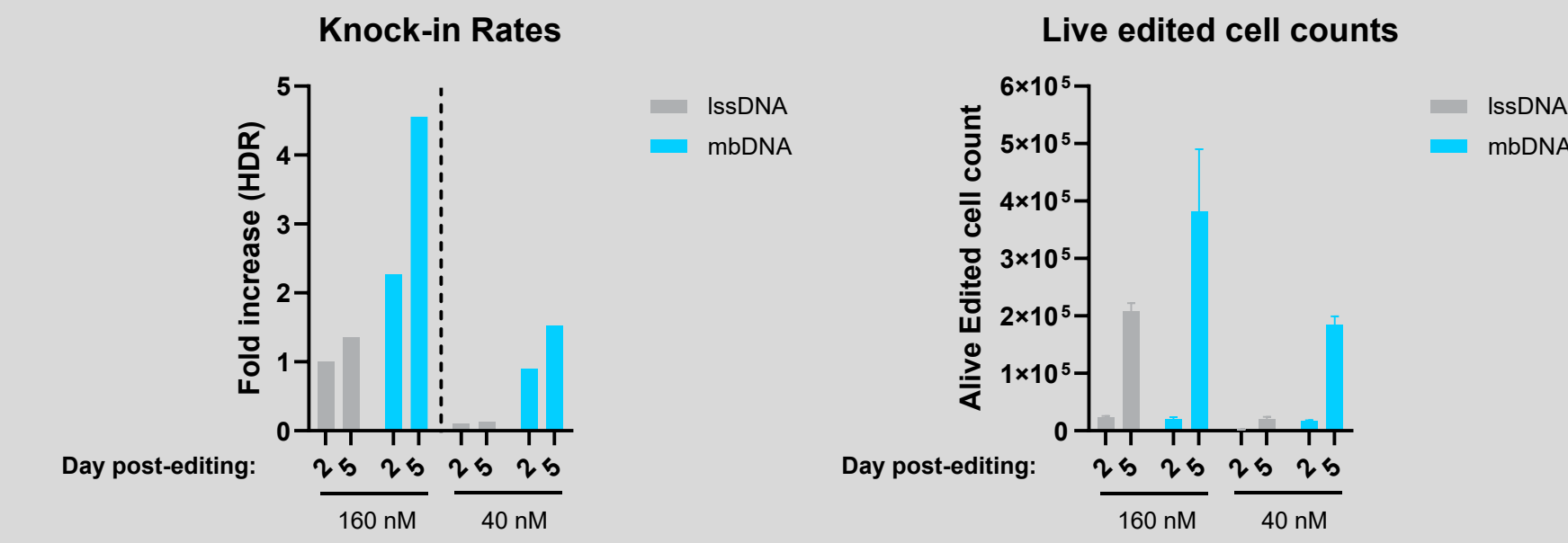

Higher KI rates and live edited cell counts

Enables efficient multiplexing

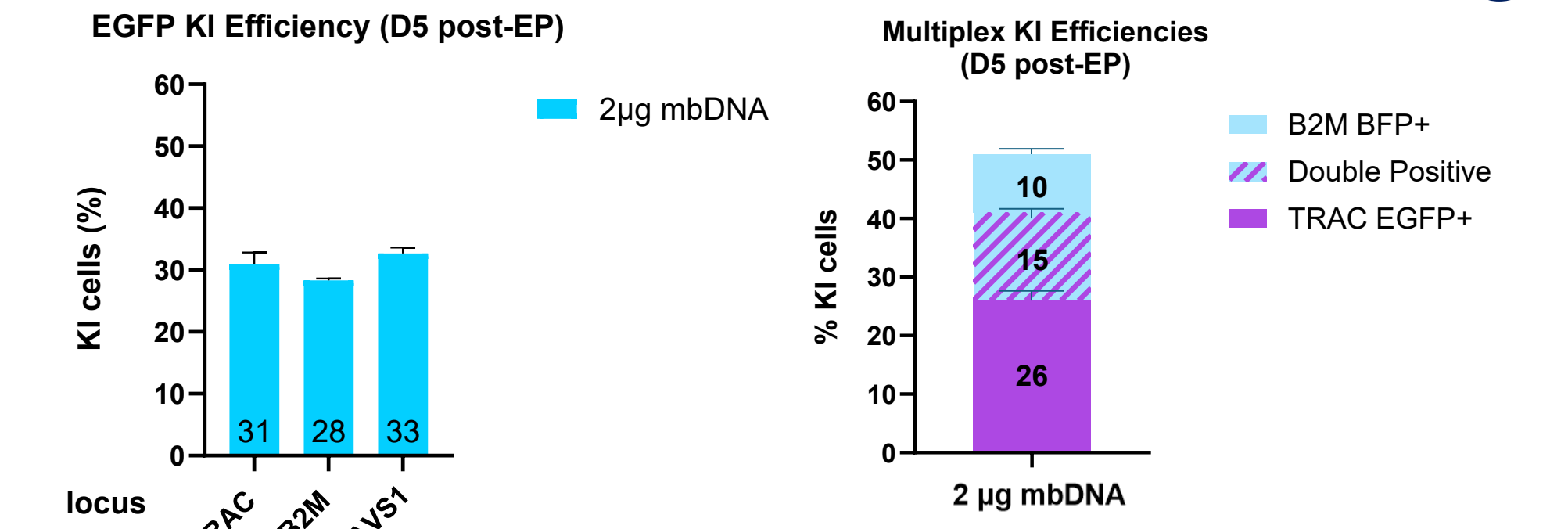


mbDNA delivers similar knock-in at the **TRAC**, **B2M** and **AAVS1** loci in T cells

Consistent editing at **multiple loci** allows efficient **multiplexing** and generation of **double-positive** T cells

mbDNA is a functional episomal expression vector

Luciferase expression in HepG2 cells

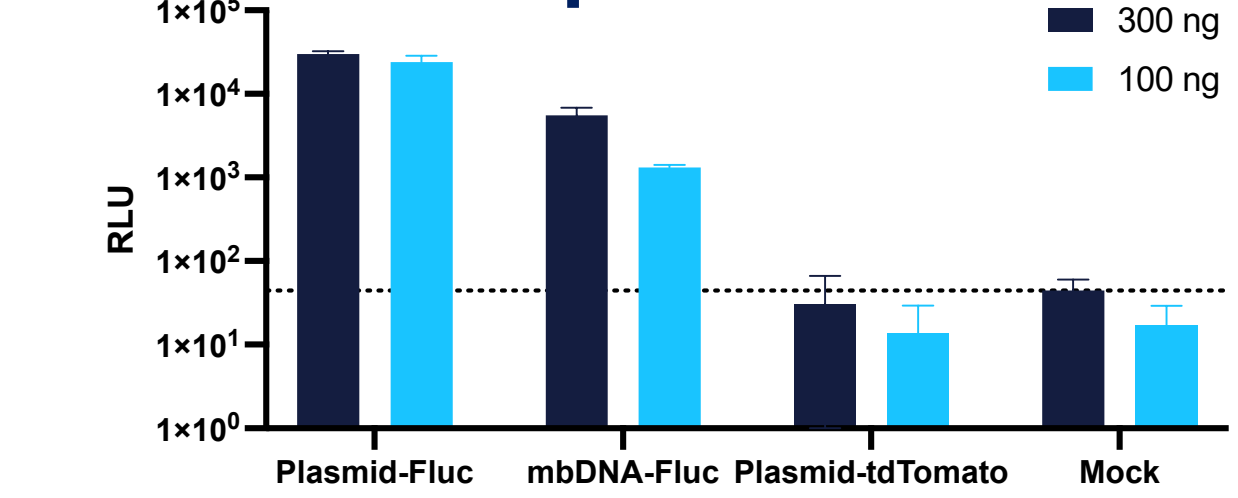


- A non-integrative mbDNA generates firefly luciferase **transgene expression** sufficiently above background in the liver cell line HepG2
- This suggests mbDNA could be used as a vector for **episomal expression** in Non-Viral Gene Therapy (NVGT), helping to **improve toxicities from double-stranded DNA** *ex vivo/in vivo*

Authors: Elizabeth Del Greco, Margarita Krivega, Julia Poniatowski and Ivan Krivega
SonoThera, Inc.

SONOTHERA™

Non-HDR Applications

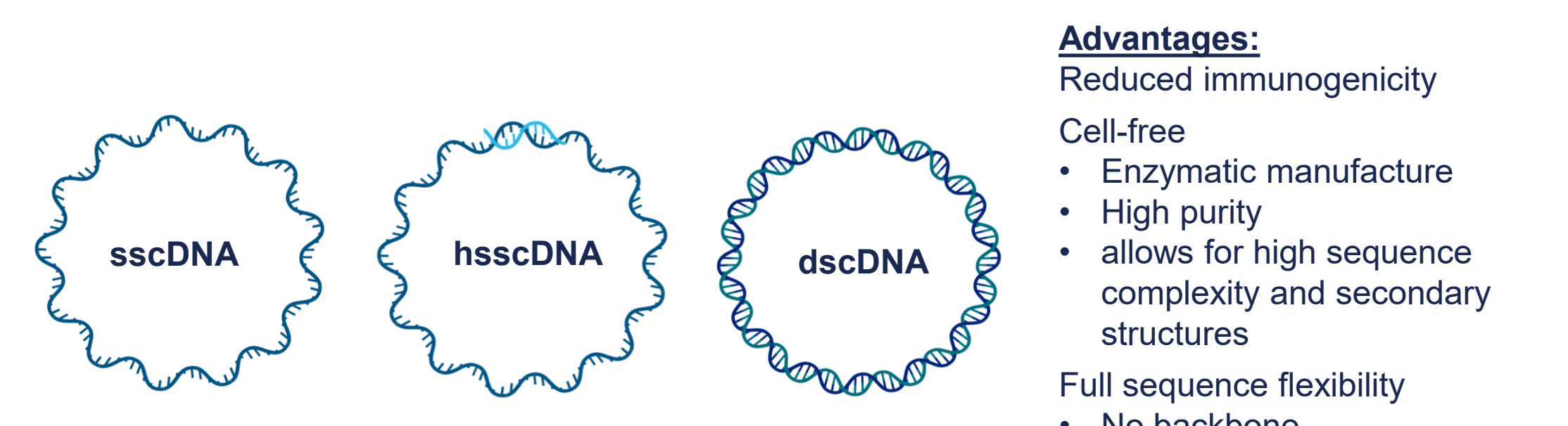

Architecture	Unique Feature	Application
Megabulb DNA (mbDNA)	Double stranded DNA stem with single stranded circle 'bulb'	Homology-directed repair (HDR), episomal expression
Double-stranded circle (dscDNA)	Fully double stranded circular molecule where every base pair is user defined	Transposases, recombinases, integrases, homology-independent targeted integration (HITI), episomal expression
Single-stranded circle (sscDNA)	Fully single stranded circular molecule where every base pair is user defined	
Hybrid single-stranded circle (hsscDNA)	User defined double stranded and single stranded regions Phosphorothioates to protect DNA ends from degradation	